

Spare parts – BioCompact II

RR 610 H




Description :

Spare parts list for BioCompact II
RR 610 H

Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
12.01.09	JP	14.10.24	Jolu	026		


Exploded view RR 610

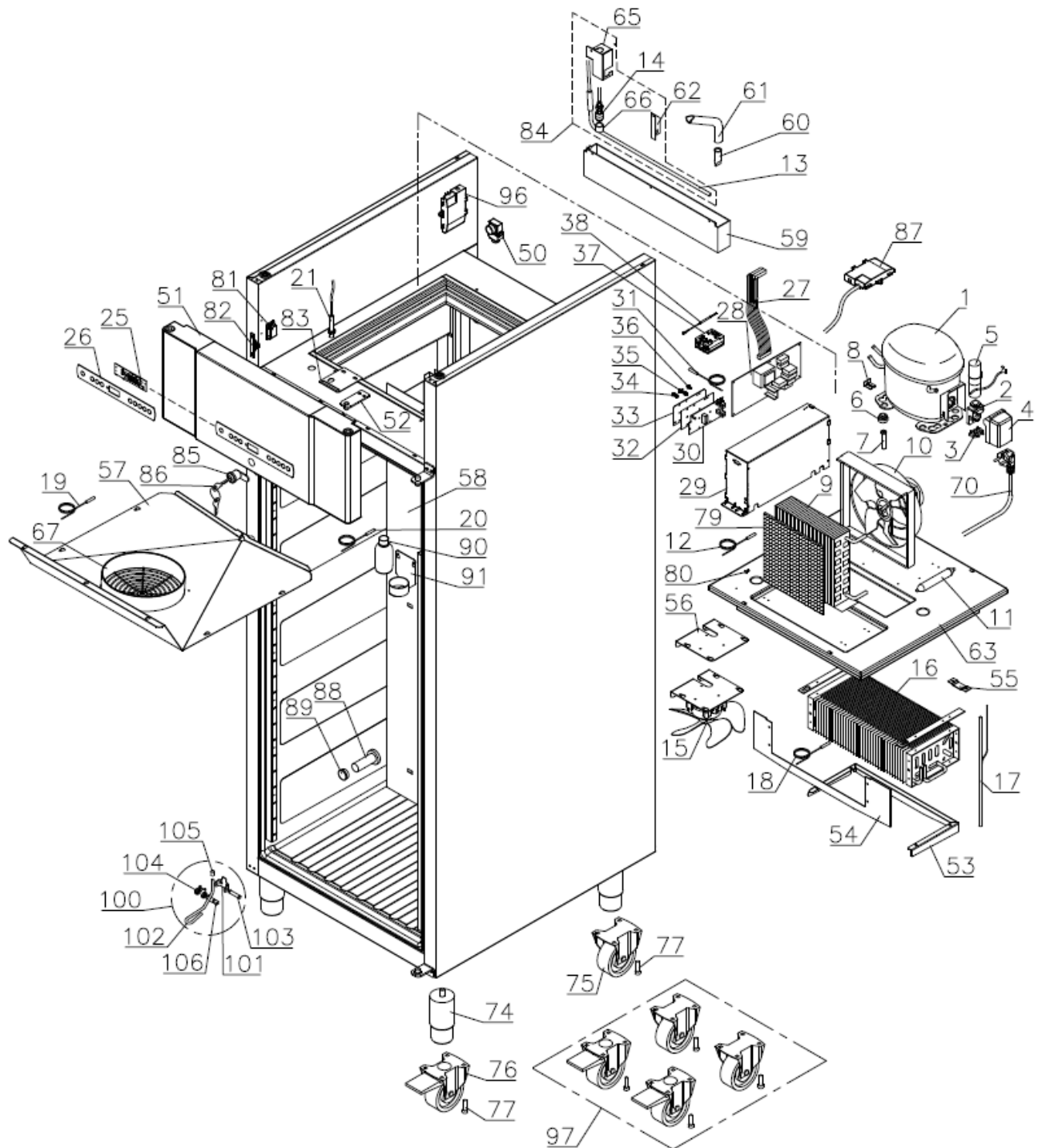

Technical information					765041295	
<u>Description :</u>						
Spare parts list for BioCompact II RR 610 H						
Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
12.01.09	JP	14.10.24	Jolu	026		


Door

0508-4314

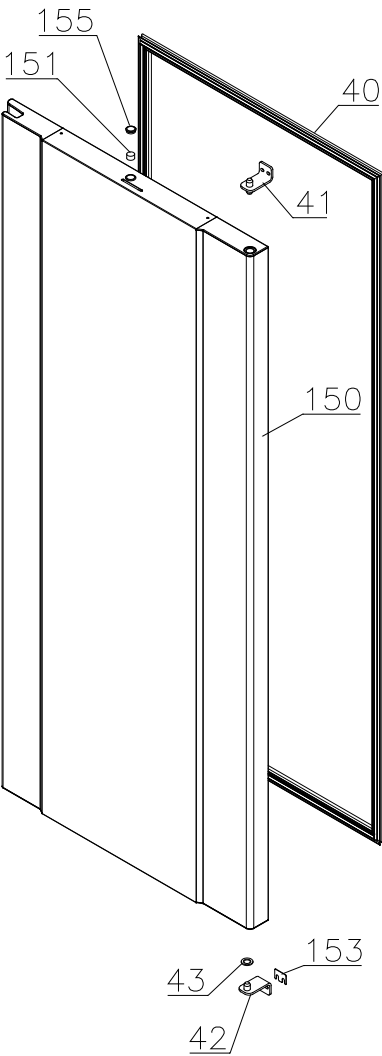

Technical information					765041295	
<u>Description :</u> Spare parts list for BioCompact II RR 610 H						
Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
12.01.09	JP	14.10.24	Jolu	026		

4314-

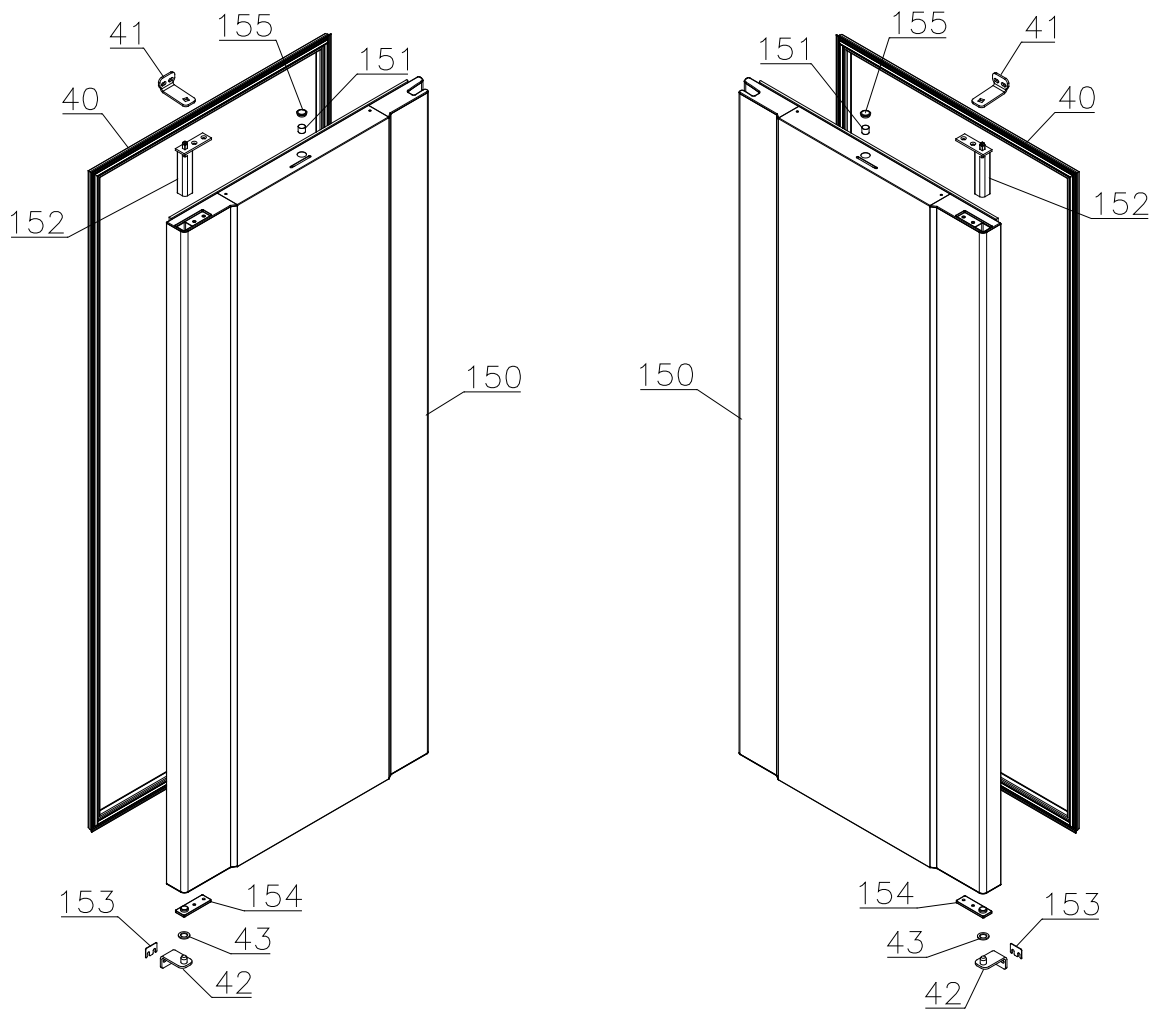


Description :

Spare parts list for BioCompact II
RR 610 H


Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
12.01.09	JP	14.10.24	Jolu	026		

Chart recorder parts

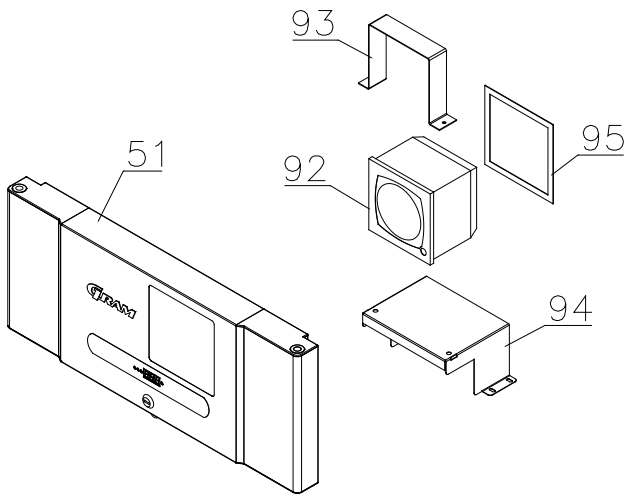

Interior parts

Shelves:

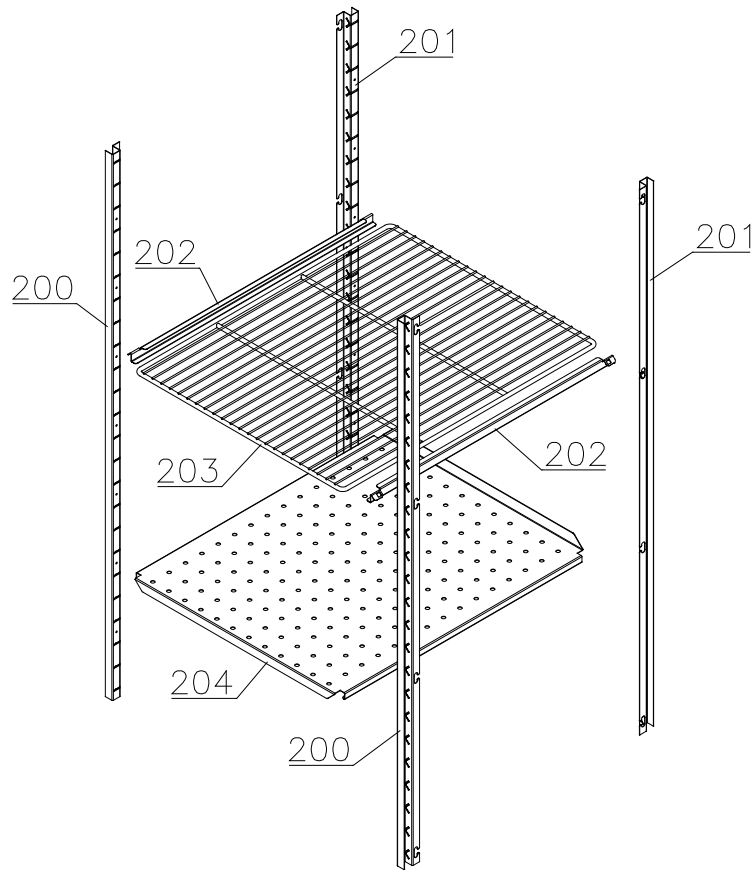


Description :

Spare parts list for BioCompact II
RR 610 H

Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
12.01.09	JP	14.10.24	Jolu	026		

Drawers:

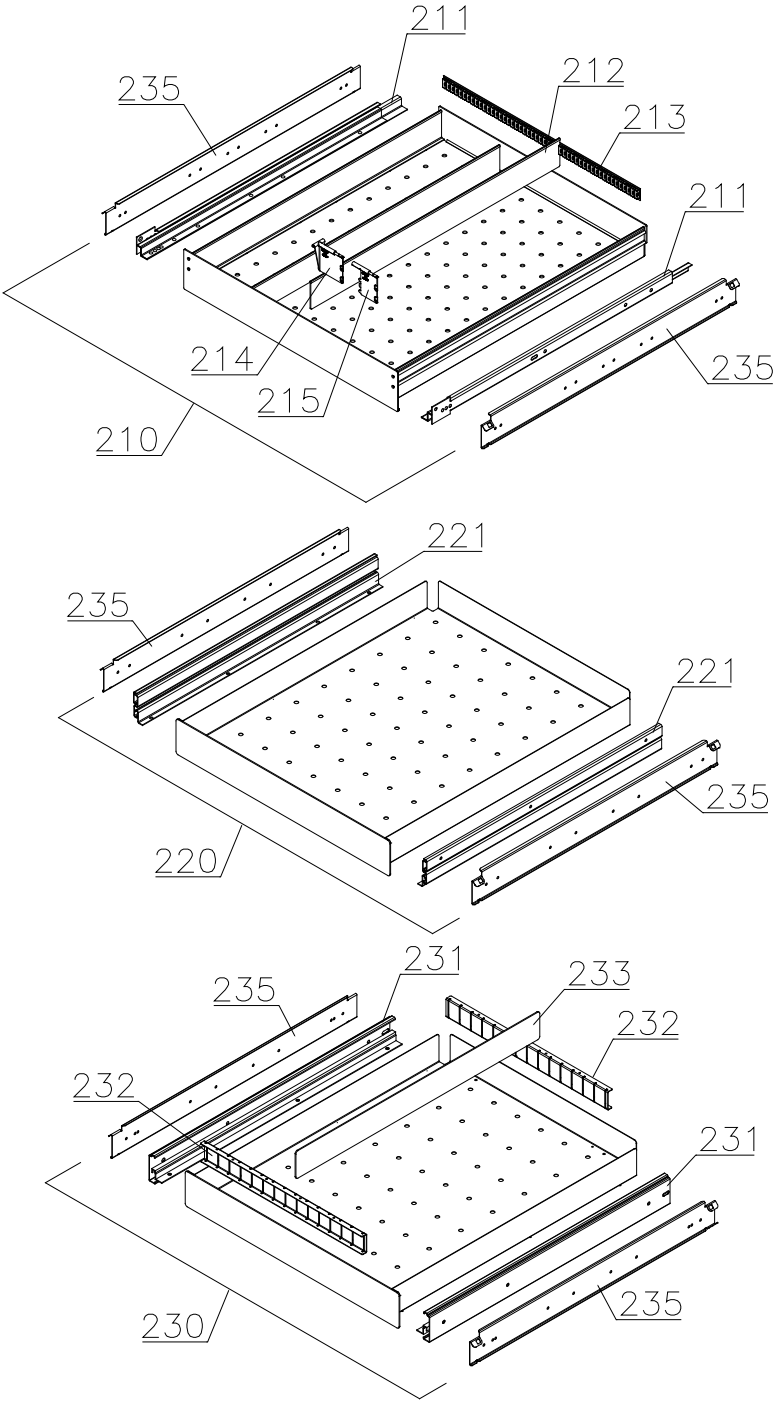

<i>Technical information</i>					765041295	
<u>Description :</u> Spare parts list for BioCompact II RR 610 H						
Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
12.01.09	JP	14.10.24	Jolu	026		

Item no.	Description	Gram part no.	week/year introduced	week/year deleted
----------	-------------	---------------	----------------------	-------------------

Essential spare parts

1	Compressor FR 7.5 G - 50/60 Hz	500550061	1508	
2	Starting relay	410589015	1508	
3	Cord relief	452440297	1508	
4	Cover	452440293	1508	
5	Start capacitor	451720018	1508	
6	Rubber foot	760080519	1508	
7	Screw for compressor	423200210	1508	
8	Fastening clip	760082367	1508	
9	Condenser	612900500	1508	
10	Condenser fan guard compl.	765080189	1508	
11	Filter drier	762900157	1508	
12	Condenser sensor	453531500	1508	
13	Re-evaporation heater	761250778	1508	
14	Float switch	765072049	1508	
15	Evaporator fan complete	761221902	1508	3921
15	Evaporator fan complete	761221904	3921	
16	Evaporator	761220332	1508	
17	Heat exchanger	761221403	1508	
18	Defrost sensor	453531501	1508	
19	Cabinet sensor	453531501	1508	
20	E-sensor	453531504	1508	
21	Door switch	450400113	1508	
25	Controller display part	612900056	1508	
26	Label with keys	765041752	1508	
27	Band cable	453582046	1508	
27	Band cable for LTP	765072958	3519	
28	Electronic controller K2+	761256120	1508	
29	Electrical box complete	761232111	1508	
30	LTP controller	453580515	3519	
31	LTP sensor	453531526	3519	
32	PCB protection sheet	765072698	0718	
33	Screening plate	765073271	0718	
34	Dual locking spacer H=7.9 mm	451119525	0718	
35	Dual locking spacer H=15.9 mm	451119526	0718	
36	Push spacer H=7.9 mm	451119527	0718	
37	Solid state relay	450200300	4717	
38	Cable to SSR incl. resistance	765071445	1023	
38	Cable to SSR incl. resistance "LTP"	765071449	1023	

Technical information					765041295	
<u>Description :</u> Spare parts list for BioCompact II RR 610 H						
Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
12.01.09	JP	14.10.24	Jolu	026		

Item no.	Description	Gram part no.	week/year introduced	week/year deleted
40	Door gasket	761222863	1508	
41	Hinge, top	761220385	1508	4314
41	Hinge, top, self-close	761222802	4314	
41	Hinge, top, self-close	761222803	4314	
42	Hinge, bottom	761220387	1508	4314
42	Hinge, bottom, self-close	761222819	4314	
43	Washer	760250524	1508	4314
43	Washer	421708003	4314	

<i>Technical information</i>					765041295	
<u>Description :</u> Spare parts list for BioCompact II RR 610 H						
Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
12.01.09	JP	14.10.24	Jolu	026		

Item no.	Description	Gram part no.	week/year introduced	week/year deleted
----------	-------------	---------------	----------------------	-------------------

Other spare parts

50	Air relief valve	765081206	1508	
51	Front panel stainless	761231403	0508	
51	Front panel white	761231404	0508	
51	Front panel f. chart recorder stainless	761231451	4400	
51	Front panel f. chart recorder white	761231452	4400	
52	Hinge for front panel	761220388	1508	
53	Evaporator cover, side/back	761220326	1508	
54	Evaporator cover, front	761220325	1508	
55	Suspension bracket	760250295	1508	
56	Fan holder	761220369	1508	
57	Drip water tray	761221430	1508	
58	Air distribution plate	761220331	1508	
59	Re-evaporation tray	765081149	1508	
60	Water trap	760082891	1508	
61	Drain pipe	765081153	1508	5017
61	Drain pipe	765081253	5017	
62	Heater bracket	761255445	1508	2714
62	Heater bracket	763250002	2714	
63	Insulation slab	761221734	1508	
64	Screw for fan blade	612900100	1508	
65	Float switch bracket	761255446	1508	2714
65	Float switch bracket	763250002	2714	
66	Plastic cap	450450080	1508	
67	Fan guard	765081239	1508	
70	Lead and plug, DK 3-pin	765072568	1508	
70	Lead and plug, Schuko	765072566	1508	
70	Lead and plug, Switzerland	765070163	1508	
70	Lead and plug, GB	765072567	1508	
74	Leg, stainless H=135-200	765080115	1508	
75	Fixed wheel, H=125	480270520	1508	
76	Castor wheel, with brake	480260050	1508	
77	Screw M8 x 12	422715000	1508	
79	Condenser filter	761114416	1508	
80	Attachment spring	762110375	1508	289
80	Attachment plug	765081246	2809	
81	Front panel close catch	761220375	1508	
82	Front panel close keeper	761253301	1508	
83	Bracket for lock	761220376	1508	
84	Re-evap. heater kit complete	763250010	1508	

<i>Technical information</i>					765041295	
<u>Description :</u> Spare parts list for BioCompact II RR 610 H						
Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
12.01.09	JP	14.10.24	Jolu	026		

Item no.	Description	Gram part no.	week/year introduced	week/year deleted	
85	Door lock	765080092	1508		
86	Key no. 1 (25010)	765080340	1508		
86	Key no. 2 (25020)	765080341	1508		
86	Key no. 3 (25030)	765080342	1508		
86	Key no. 4 (25040)	765080343	1508		
86	Key no. 5 (25050)	765080344	1508		
86	Master key	765080385	1508		
87	Box for alarm connection	761233402	1508		
88	Insulation plug	765080563	1508		
89	Plastic cap	501020023	1508		
90	Liquid container	765080384	1508		
91	Holder for liquid container	761256497	1508		
92	Chart recorder	761256438	1508		
93	Bracket for chart recorder	761256439	1508		
94	Base for chart recorder	761231447	1508		
95	Cover for chart recorder	761236452	1508		
96	Electrics cover	765070143	1508		
97	Set of castor wheels	818899905	1508		
100	Foot pedal complete	819581053	0911		
101	Bracket for pedal	761224010	0911		
102	Pedal arm	761224002	0911		
103	Spindle	761224007	0911		
104	Spring	761224005	0911		
105	Plastic cap	581400021	0911		
106	Bush	761224008	0911		
150	Door complete	stainless	761230917	0508	4314
150	Door complete	white	761230916	0508	4314
150	Door complete, self-close, right	stainless	761232903	4314	4024
150	Door complete, self-close, right	stainless	762234014	4124	
150	Door complete, self-close, left	stainless	761232913	4314	4024
150	Door complete, self-close, left	stainless	762234015	4124	
150	Door complete, self-close, right	white	761232904	4314	4024
150	Door complete, self-close, right	white	762234012	4124	
150	Door complete, self-close, left	white	761232914	4314	4024
150	Door complete, self-close, left	white	762234013	4124	
151	Magnet for door switch		450401002	1508	
152	Self-close device		763251745	4314	
153	Spacer for hinge		761258621	0508	
154	Hinge bearing complete		763251755	4314	
155	Plastic cap, white		760082756	0508	
155	Plastic cap, grey		765081220	0508	

<i>Technical information</i>					765041295	
<u>Description :</u> Spare parts list for BioCompact II RR 610 H						
Compiled.		Latest revision.			Deleted.	
Date	Name	Date	Name	Rev.no	Date	Name
12.01.09	JP	14.10.24	Jolu	026		

Item no.	Description	Gram part no.	week/year introduced	week/year deleted
----------	-------------	---------------	----------------------	-------------------

Interior equipment:

200	Wall rail, front	761220335	1508	
201	Wall rail, back	761220336	1508	
202	Shelf support rail	761256455	1508	
203	Grey wire shelf	818880008	1508	
204	Perforated shelf, stainless	761230083	1508	
210	Alu drawer, complete	761230362	2210	
211	Set of sliding rails	761234055	2210	
212	Divider rail	761230016	2210	
213	Holder for divider rail	761230063	2210	
214	Plastic divider 1	761230021	2210	
215	Plastic divider 2	761230022	2210	
220	Stainless drawer 30 kg, complete	761231232	1508	
221	Set of sliding rails	761234000	1508	
230	Stainless drawer 40 kg, complete	761231202	1508	
231	Set of sliding rails	761234058	1508	
232	Divider rail	761230305	1508	
233	Divider	761230308	1508	
235	Support for sliding rail	761230201	1508	3310
235	Support for sliding rail, right	761230278	3310	
235	Support for sliding rail, left	761230281	3310	